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## ⚠ Warning

BEFORE USING MYPALLETIZER READ ALL INSTRUCTIONS AND CAUTIONARY MARKINGS IN THIS MANUAL

1. Do not expose the product to rain or moisture to reduce fire or shock hazard.
2. Do not place the product in or near fire.
3. Do not leave the product in a car in hot or humid weather.
4. Do not disassemble, crush or pierce the product.
5. Do not expose the product to excessive shock such as dropping from a high place.
6. Do not expose the product to high temperatures above 60 °C (140 °F).

## ⚠ Attention

⚠ Regarding the operation and secondary development of myPalletizer, please read and download the related files before using it.

1、Gitbook 2、User Manual 3、Development Manual

Official Website: <https://www.elephantrobotics.com/en/mypalletizer-en1/>

## Zero Basic Entry Level-Desktop Robotic Arm

The new series of myPalletizer elephant palletizing robot arm, which is jointly produced with M5STACK, is a fully packaged lightweight four-axis palletizing robot arm. The overall finless design is compact and easy to carry.

myPalletizer body weight 960g, load 250g, working radius 260mm; specially designed for makers and education, easy to use, rich expansion interfaces, unlimited development possibilities; a variety of AI suit options, machine vision, composite robots, etc., learn about the cutting-edge technology of learning robots Knowledge inspires innovative thinking.

### The fin design, the appearance is fully wrapped

The optimal space-removing fin design concept that can be loaded into a backpack subverts the traditional link-type educational four-axis robotic arm.


### LED dual display

The robotic arm carries two display screens, which is easy to expand the interactive output of the application, and the gameplay is unlimited.

### LEGO ecology, compatible with all my series accessories

The patented Lego hole design is shared globally, and the my series hardware ecological platform concept is implemented, and the end accessories are plug and-play.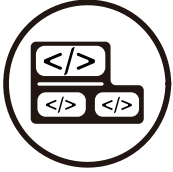

### Zero foundation, easy to get started


It is easy to get started and use, and the graphical programming language allows you to easily start the journey of using the robotic arm.

### Support ROS ecology

It is developed using the global mainstream robot communication framework ROS, and supports simulation, control and algorithm verification in a virtual environment, which reduces the requirements for the experimental environment


## myPalletizer - Size and Working Range Diagram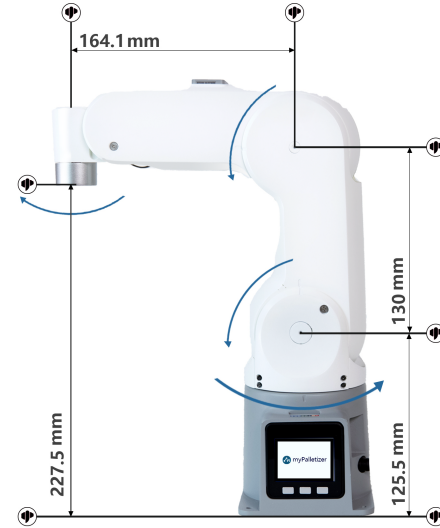

## Parameter

Name	myPalletizer
Model	myPalletizer 260 M5
Degree of Freedom	4
Repeatability	± 2mm
Payload	250g
Weight	960g
Working radius	260mm
Material	Photosensitive resin SLA
Power Input	8 ~ 12V 5A
Motor type	High precision magnetic encoder servo motor
Max speed	120°/s
Control	M5Stack Basic
Communication	USB / type-C


## M5STACK Control Board Pin Map

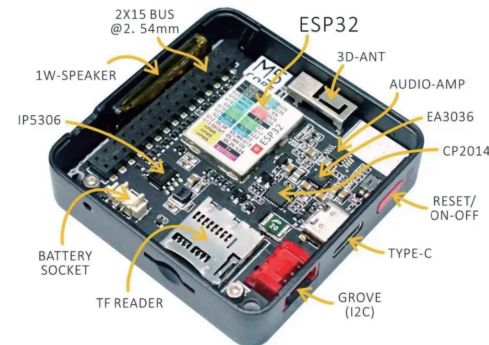
M5 Basic side pin diagram1

Atom diagram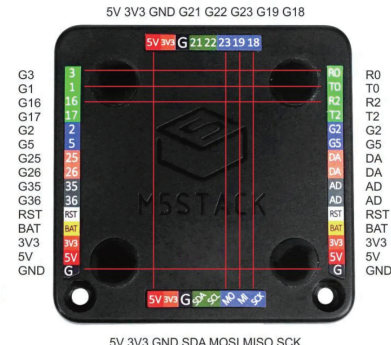


M5 Basic side pin diagram1

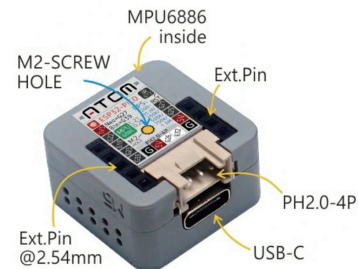Atom pin diagram

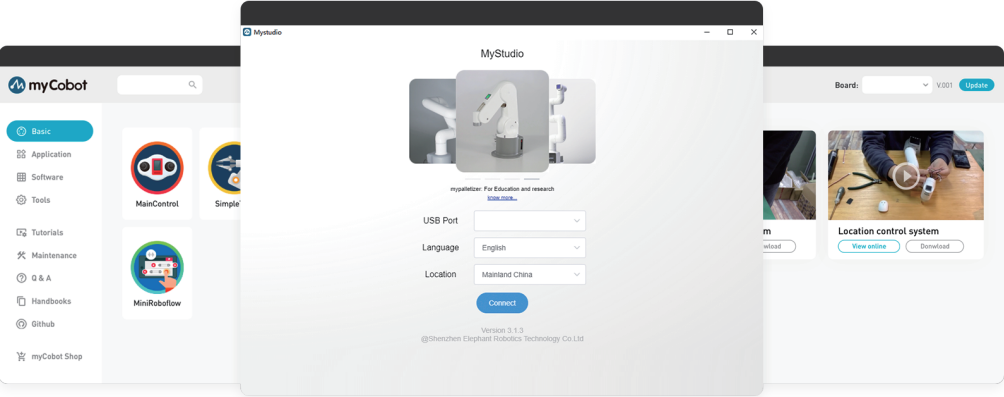myStudio is a one-stop platform for robots


myStudio integrates my series software and various materials. The main functions of myStudio are: 1) Update the firmware; 2) Provide video tutorials on how to use the robot; 3) Provide maintenance and repair information (such as video tutorials, Q&A, etc.).

Please download the latest version of myStudio to use.

The download link is as follows:  
Officail website: <https://www.elephantrobotics.com/en/downloads/>  
Github: <https://github.com/elephantrobotics/MyStudio/>

Burn Table

Please use myStudio software to burn firmware. Development environments that support the secondary development of myPalletizer are: myBlockly, RoboFlow, Arduino, ROS, python, etc.

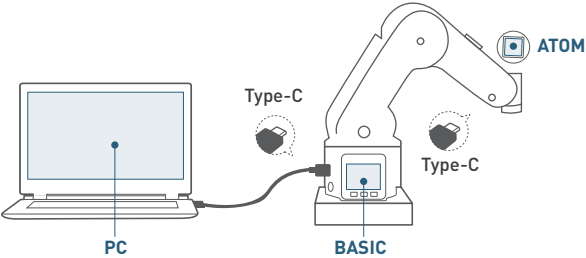Development Environment	Library on PC	Basic Firmware	Atom Firmware
Default Program	N/a	mainControl	atomMain
Visual Programming	myBlockly	Transponder	atomMain
RoboFlow Industrial Programming Software	RoboFlow	Transponder	atomMain
Arduino Maker!	Arduino IDE + M5Stack Lib + MycobotBasic Lib	All Exapmles	atomMain
API on Desktop	Python/ C+	Transponder	atomMain
ROS Development	ROS	Transponder	atomMain
USB/TxRx0(G1/G3)	Read Protocol	Transponder	atomMain
BlueTooth	Read Protocol	BT_Transponder	atomMain
phoneApp	Mobile phone Android/iPhone	BT_Transponder	atomMain

# myPalletizer Accessory

Adaptive Gripper

Camera Flange

Suction Pump

G Base

Elephant Robotics are targeted at robotic collaboration applications, making “my-series” product line. For new information about the accessories, Follow us on Shopify and Twitter.  
**Shopify:** <https://shop.elephantrobotics.com/>  
**Twitter:** @cobotMy

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Purchaser \_\_\_\_\_ Order No. \_\_\_\_\_ Phone \_\_\_\_\_

Address \_\_\_\_\_ Logistics Receipt Date \_\_\_\_\_

Product problem description(Required):

\_\_\_\_\_

If you need to apply for warranty service, please contact our customer service to confirm the detailed information. After confirmation, please fill in the card and send it back together with the product and the attached invoice. **Note: Our company reserves the right to explain and modify the warranty card of this product within the scope of the law.**

- Return service is limited to goods not opened within 7 days after the receipt date of logistics of the products. The freight or other risks incurred in return shall be borne by the customer.
- Customers should provide the purchasing invoice and warranty card as the warranty certification when a warranty is being asked.
- Elephant Robotics will be responsible for the hardware faults of products caused by the normal using during the warranty period.
- The warranty period starts from the date of purchase or the receipt date of the logistics.
- The faulty parts from the products will be owned by Elephant Robotics, and the appropriate cost will be charged if necessary.

If you need to apply for warranty service, please contact our customer service first to confirm the detailed information.

servo motor	
Warranty Period	Warranty Services
≤1 months	Elephant Robotics offers a free new servo motor and bear the freigh.
1-3 months	Elephant Robotics offers a free new servo motor, customs shall bear the freight.
≥3 months	Customers need to buy it themselves.
Electrical Parts (M5 Hardware)	
≤3 months	Customers need to send it back after disassembly, Elephant Robotics shall send a new one for free and bear the freight out and home.
3-6 months	Customers need to send it back after disassembly and bear the freight out and home, Elephant Robotics shall send a new one for free.
≥6 months	Customers need to buy it themselves.
Structure Parts, including Shell Parts	
≤1 year	Elephant Robotics offers free new components once, customs shall bear the freight.
≥1 year	Customers need to buy it themselves.

During the warranty period of the delivered product, the company only repairs the malfunctions that occur during normal use of the robot for free. However, in the following cases, the customer will be charged for repairs (even during the warranty period):

- Damage or malfunction caused by incorrect use and improper use different from the contents of the manual.
- Failure caused by unauthorized disassembly by the customer.
- Damage caused by improper adjustment or unauthorized repairs.
- Damage caused by natural disasters such as earthquakes and floods.

Therefore, please strictly follow the instructions in this manual and related manual to operate the robot.

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© 2020 Shenzhen Elephant Robotics Technology Co.Ltd