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### ⚠ Warning

BEFORE USING MYPALLETIZER READ ALL INSTRUCTIONS AND CAUTIONARY MARKINGS IN THIS MANUAL

1. Do not expose the product to rain or moisture to reduce fire or shock hazard.
2. Do not place the product in or near fire.
3. Do not leave the product in a car in hot or humid weather.
4. Do not disassemble, crush or pierce the product.
5. Do not expose the product to excessive shock such as dropping from a high place.
6. Do not expose the product to high temperatures above 60 °C (140 °F).

### ⚠ Attention

⚠ Regarding the operation and secondary development of ultraarm-p340, please read and download the related files before using it.

1、Gitbook 2、User Manual 3、Development Manual

Official Website: <https://www.elephantrobotics.com/en/ultraarm-p340-en/>

## ultraArm P340

### For Robot Starters - multifunctional desktop robot arm

ultraArm is a small desktop robotic arm. It is designed with classic metal structure and occupies only an area of A5 paper. It is equipped with high-performance stepper motor, and it owns  $\pm 0.1\text{mm}$  repeated positioning accuracy and high stability. UltraArm can achieve the precise writing and drawing, laser engraving, and can be freely matched with multiple accessories such as slide rail, conveyor belt, myAGV and cameras. It can complete different practical training projects according to different needs, and is widely used in many fields such as higher education, scientific research, vocational application education, etc.

## Vision education kit

These kits compose of open source AI algorithm, support related teaching application scenarios, identifying and sorting objects, on slide rail or conveyor belt. There are a variety of sets with static vision or dynamic vision in a one-stop platform to help users get started quickly.

## Multiple programming environments supporting


Drag-and-drop programming visualization software: myBlockly; Provides the interface for Python developing; Provides open-source code to help users deeply learn robotics.


OpenCV


Python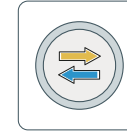


Inverse kinematics

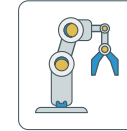Principles of  
Robot Control

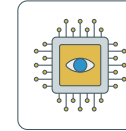Visual recognition

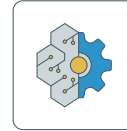End-effector use


## Hardcore configuration

It is equipped with stepper motor, and it owns  $\pm 0.1\text{mm}$  repeated positioning accuracy; It supports 24/7 highly intensive operation; It is compact and tough with classic metal structure.

## DIY kit

This is a one-click track generation in visualization software with operation plate, pen clip, laser engraving head. Users will enjoy warm moments of online production, nameplate and phone case DIY production, AI drawing.

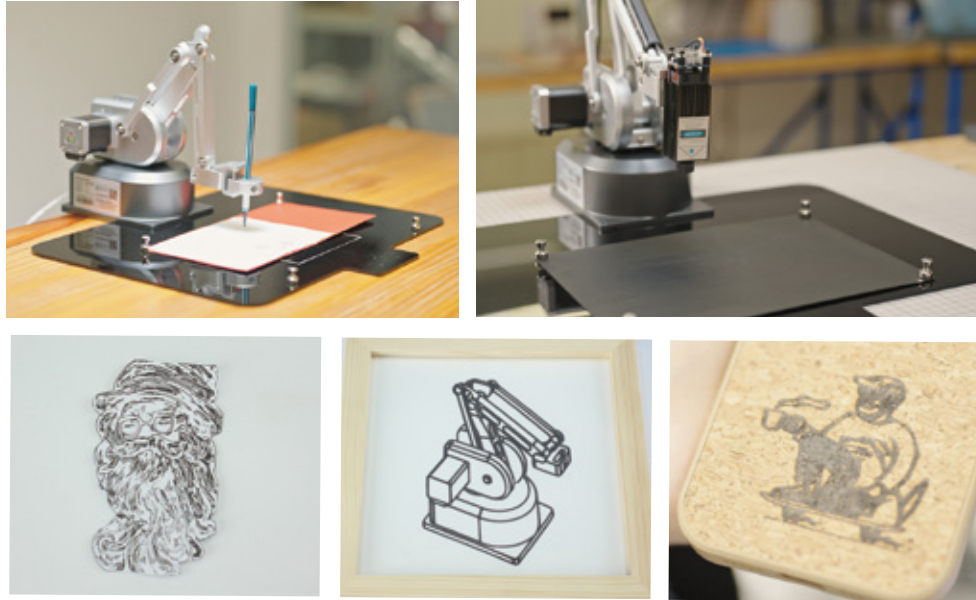## ultraArm P340 Size and working space

## Parameter

Name	ultraArm P340	Adapters	100-240VAC 50-60Hz
Degree of freedom	3DOF / 4DOF	Interfaces	Laser engraving interface
Payload	650g		Adaptive gripper interface
Working radius	340mm		Suction pump interface
Repeatability	±0.1mm		Burning switch
Weight	2.9kg	TCP terminal	100 mm/s
Working lifespan	5000h, support 7x24h	Installation mode	Desktop level
		Power module	High performance stepper motor
Power input	12V 5A	Elephant Luban,	
Available angle of joints	J1 -150° ~ +170°	Programming mode	Graphical programming myBlockly,
	J2 -20° ~ 90°		Python, ROS2
	J3 -5° ~ 75°	Temperature	0°~45°
Noise	60 dB	Ambient humidity	5%-80%
Working speed	100mm/s		

## Introduction of interfaces

Title	Sign-Name	Function	Remark
Burn	Key Switch	Burning function is on/off	When the burning function is turned on, the serial port connection will restart the development board. If you do not need to restart, please turn off the burning function
Laser	PWM GND 12V	Laser engraving interface	PWM controls laser engraving power
Gripper	DATA 8.4V GND	Gripper interface	DATA control claw/servo data
Switch	OUT1 OUT2 5V GND	Suction pump interface	5V-OUT-PIN output: control solenoid valve and suction pump switch

ultraArm P340 Accessory

Suction pump

Adaptive gripper

Laser engraving head

Servo

Pen holder

Lego connectors

Elephant Robotics are targeted at robotic collaboration applications, making “ultra-series” product line. For new information about the accessories, Follow us on Shopify and Twitter.  
**Shopify:** <https://shop.elephantrobotics.com/>  
**Twitter:** @cobotMy

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Purchaser \_\_\_\_\_ Order No. \_\_\_\_\_ Phone \_\_\_\_\_

Address \_\_\_\_\_ Logistics Receipt Date \_\_\_\_\_

Product problem description(Required):

\_\_\_\_\_

If you need to apply for warranty service, please contact our customer service to confirm the detailed information. After confirmation, please fill in the card and send it back together with the product and the attached invoice. **Note: Our company reserves the right to explain and modify the warranty card of this product within the scope of the law.**

- Return service is limited to goods not opened within 7 days after the receipt date of logistics of the products. The freight or other risks incurred in return shall be borne by the customer.
- Customers should provide the purchasing invoice and warranty card as the warranty certification when a warranty is being asked.
- Elephant Robotics will be responsible for the hardware faults of products caused by the normal using during the warranty period.
- The warranty period starts from the date of purchase or the receipt date of the logistics.
- The faulty parts from the products will be owned by Elephant Robotics, and the appropriate cost will be charged if necessary.

If you need to apply for warranty service, please contact our customer service first to confirm the detailed information.

Warranty policy	
≤1 year	Elephant Robotics offers free new components once, customs shall bear the freight.
>1 year	Customers need to buy it themselves.

During the warranty period of the delivered product, the company only repairs the malfunctions that occur during normal use of the robot for free. However, in the following cases, the customer will be charged for repairs (even during the warranty period):

- Damage or malfunction caused by incorrect use and improper use different from the contents of the manual.
- Failure caused by unauthorized disassembly by the customer.
- Damage caused by improper adjustment or unauthorized repairs.
- Damage caused by natural disasters such as earthquakes and floods.

Therefore, please strictly follow the instructions in this manual and related manual to operate the robot.